



New UI FAQ #15

How do I Result a Manual Test?

To Result a Manual Test:

1. Click on **Result** within the Universal Add button or click on the **Add** icon in the Results data table:

Results / Flagged

Status: 2 Sample Type: Patient Result Time: Previous

STATUS	RESULT TIME	DEVICE TYPE	SAMPLE ID	OPERATOR	PATIENT	LOCATION	INSTITUTION	FLAGS	DEVICE	OUTBOUND FAI
Evaluate	1/12/2022 9:20:00 PM	COVID-19	58339	COVID OP (01220857...		2 Tower (Clifton Avenue...	Clifton Avenue Medica...	Unknown Test Material		
Evaluate	1/12/2022 2:30:00 PM	Pregnancy	951357	COVID OP (01220857...		2 Tower (Clifton Avenue...	Clifton Avenue Medica...	Uncertified Operator (...)		
Evaluate	1/12/2022 1:03:00 PM	COVID-19	572055	COVID OP (01220857...		2C (Clifton Avenue Me...	Clifton Avenue Medica...	Unknown Test Material		
Evaluate	1/12/2022 12:17:00 PM	COVID-19	581663	COVID OP (01220857...		3E (Clifton Avenue Me...	Clifton Avenue Medica...	Unknown Test Material		
Evaluate	1/12/2022 11:37:00 AM	Pregnancy	3745	Robin Vu (82815)		ER (Clifton Avenue Me...	Clifton Avenue Medica...	Expired Test Material		
Evaluate	1/7/2022 11:35:00 AM	Pregnancy	5294	Robin Vu (82815)		2C (Clifton Avenue Me...	Clifton Avenue Medica...	Expired Test Material		
Evaluate	1/7/2022 10:44:00 AM	Fecal Occult	1399	Robin Vu (82815)		2E (Clifton Avenue Me...	Clifton Avenue Medica...	Uncertified Operator (...)		

Add Result

2. Complete the **General** section, which includes selecting the **Location** the test was completed in, the **Manual Test** being resultated, and select the **Sample Type**, then click **Next**

Results / Add Result

1 General

Location *

2C (Clifton Avenue Medical Unit)

Manual Test *

BinaxNOW COVID-19 Ag (COVID-19)

Sample Type *

Patient

NEXT



How do I Result a Manual Test?

3. Complete the **Sample** section, filling in the required and optional fields as configured within RALS. Required fields will have an asterisk next to the title. Configuration of the manual test will vary, be sure to follow internal procedures when resulting. After completing the **Sample** details, click **Next**

2 Sample

Sample ID *

2562

No patient found
Kit Lot *

1204 (Kit)

Collection Time

Result Time *

1/13/2022 8:43 AM

Specimen Type *

Swab (COVID-19)

Q Specimen Site

BACK

NEXT

4. Enter the manual test result/s, then click **Next**

3 Values

PARAMETER	VALUE	UNIT OF MI
Control Line	Valid	
SARS-CoV-2	Negative	
Test Kit	345	

BACK

NEXT



New UI FAQ #15

How do I Result a Manual Test?

5. Comments can be added by clicking the **Add** icon. Select from the drop-down list or enter a free-text comment.

4 Comments



Free Text Comment

BACK

SAVE

6. To complete the test, click **Save**